


Installation Guide Stylet+ software and drivers for Giga-8d

1. The *Stylet+_acq G8d (styletaq v01.30.exe)* file should be downloaded and opened on your computer
<https://epgsystems.eu/downloads-install-files-manuals/category/3-installation>

NOTE: *your organization may block downloading .exe files from the internet. Ask your system administrator for permission to do so.*

2. Click on 'continue' messages, accept copy right, etc. and next, subtract the FTDI CMD Drivers. You will end up in the demo screen for data acquisition. As indicated in the bottom right corner by 'Demo device'.
3. You may either close or continue this data acquisition demo session (see Giga-8d manual).
4. Your *C:\Stylet+* folder should now at least contain the *Stylet+a.exe*, *Stylet+d.exe* and *unins000.exe* files. However, the *Stylet+d.exe* is not compatible with the Giga-8d. The proper file for Giga-8d data acquisition is *Stylet+d_V01.30.exe*.
5. Download this *Stylet+d_V01.30.exe* file now, put it in the *C:\Stylet+* folder and make a desktop shortcut of it.
6. Delete the *Stylet+d.exe* file.

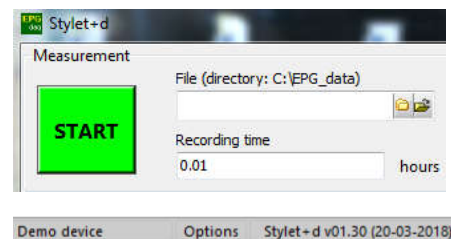

For **Giga-8d operation** (EPG recording) download and read *Manual-Giga-8d 17.pdf*

Note: In case of any installation or other problems, please contact info@epgsystems.eu