



April 19, 2017

EPG systems

WAGENINGEN, THE NETHERLANDS

-- Installation Guide for older Giga (-4 or-8) devices --

1. Downloading

Download the following 4 files on this page, and store them in C:\Stylet+\:

active.exe
add.Di710
145usbdriver.exe
Stylet setup_100516.exe

2. AD Driver Installations

1. Do not connect the USB cable of the Giga or separate Dataq device.
The Dataq AD (analog to digital) device can be inside the Giga or can be a separate device, cable connected to the Giga).
2. Go to C:\Stylet+ and execute the add.vDi710.exe (this will cause the activex.exe file to be executed as well). Follow the instructions you are prompted to choose USB interface Windaq add-ons [OK] Do not install help files. After finish select 'This program is installed correct'.
3. Then execute *145usbdriver.exe*
In addition to the AD device drivers, also the commercial Dataq acquisition software 'WindaqLite' (free version) will be installed but you will not need this. Stylet+ contains all data acquisition and analysis software needed.
After finish select 'This program is installed correct'.

3. Stylet+ EPG acquisition and analysis software installation

1. Download and run Stylet setup_100516.exe
It is easy to create desktop or toolbar shortcuts.
2. This installs all Stylet+ files in folder C:\Stylet+, creates the C:\EPG_data folder, and opens the Stylet+d recording (data acquisition) program.
3. You should see (at R-bottom) the most recent file version:
4. Close this file and run Stylet+a.exe, most recent version is:

Stylet+d v01.28 (04-05-2016)

Stylet+a v01.30 (13-04-2016)

4. Updating Stylet+ (IMPORTANT!)

Regularly check the website for more recent Stylet setup_XXXXXX.exe (XXXXXX=date) versions. Some bugs may be repaired.

* Contact your IT people if Firewall settings in institutes may prevent downloading

** Always run all installation files as 'administrator'

In case of any installation or other problems, please contact by e-mail